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Ken Girvin, Senior Optical Engineer, Photonics R&D Pingsheng Tang, PhD, Senior Scientist, Photonics R&D Terry Stange, PhD, Vice President, R&D Americas

Matthew Smith, PhD, Life Sciences Vertical Marketing Manager

- i. Since the original testing and laser selection was completed, the Hach Ultra laser vendor now claims an MTTF of more than 1,000,000 hours.
- ii. Joel C. Johnson, "Method and Apparatus for Suppressing Stray Light in Particle Detectors", US Patent 6,414,754, July, 2002.
- iii. Kenneth L. Girvin and Adam J. Reed, "Method and Apparatus for Operating a Laser in an Extinction- Type Optical Particle Detector", US Patent 7,002,682, Feb., 2006.
- iv. Akira Arimoto, et al., "Optimum Conditions for the High Frequency Noise Reduction Method in Optical Videodisc Players," published in Applied Optics, vol. 25, No. 9, May 1, 1986.
- v. Masahiro Ojima, et al., "Diode Laser Noise at Video Frequencies in Optical Videodisc Players", published in Applied Optics, vol. 25, No. 9, May 1, 1986.
- vi. K. Petermann, book entitled Laser Diode Modulation and Noise, Copyright 1988 by Kluwer Academic Publishers, Chapter 4, pp. 100 to 105.
- vii. "Long Lasting Lasers: Factors that Influence Laser Life", Particle Measuring Systems, Inc., 2006.

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